


# Photometrics® Quick Installation Guide for USB

## Software Installation


### Install the Camera-Interface Software (PVCAM®)

1. Power on your computer and insert the Photometrics flash drive into a USB port. You must log on with Administrator rights.
2. Install PVCAM.
3. Begin the automated software installation sequence.
4. Restart your computer and complete the automated software installation sequence.
5. Connect the camera to the PC.


CoolSNAP™ MYO

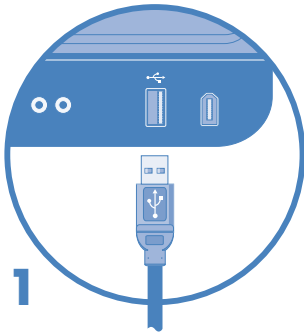


Install the USB3 PCIe card supplied with the camera.

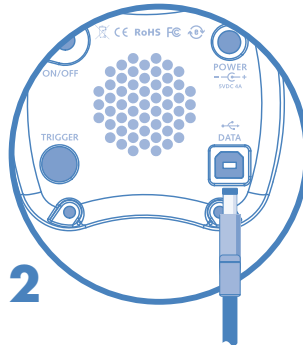

Connect the data cable to the DATA connector on the camera head. Connect the data cable to a USB port on installed card. Do not connect the cable to any other port.

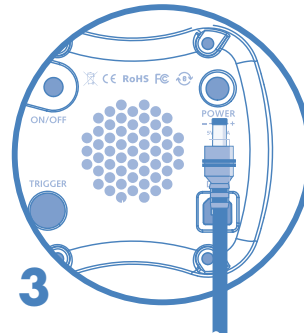

Connect the power supply to the camera's POWER connector and plug the other end of the power supply in an electrical wall outlet.

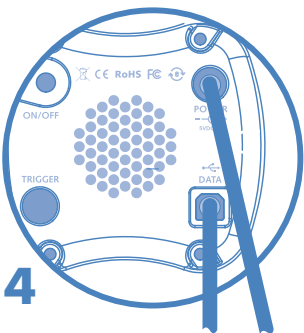

Power on your computer. Once the computer has booted up then power on the camera pushing in the ON/OFF button.

More detailed installation instructions can be found in the camera manuals on the Photometrics flash drive.

Contact Photometrics Customer Service for additional information:

+1 800.874.9789

[www.photometrics.com](http://www.photometrics.com)



Photometrics is a registered trademark.

PVCAM is a registered trademark and CoolSNAP is a trademark of Photometrics.

58-520-001 Rev A04