

Photometrics® Prime™ Software Installation Guide

Supported Operating Systems: Windows 7/8/10

Photometrics Prime Imaging Software Installation Procedure

1. Insert the included USB drive into your computer.
2. Install PVCAM 3.1.9.1. The installer is located in the root directory of the USB drive.
3. Navigate to Imaging Software > Micro-Manager
4. Install the desired version of the Micro-Manager software.
The 64-bit version is recommended.
5. Launch Micro-Manager.
6. Load the included Micro-Manager configuration file to access the camera.
The file is located in **Imaging Software > Photometrics Configuration File > PM Cam.cfg**

For a Prime sCMOS camera, use the Prime Configuration File located in

Imaging Software > Prime Configuration File > Prime.cfg

- For further system control, the configuration file will need to be updated through the Micro-Manager Hardware Configuration Wizard.

Contact Photometrics Customer Service
for additional information:

+1 800.874.9789

www.photometrics.com



Photometrics and Roper Scientific are registered trademarks of Roper Scientific, Inc. Prime and PVCAM are trademarks of Roper Scientific, Inc.

Windows is a registered trademark of Microsoft Corporation in the United States and other countries. Other brand and product names are the trademarks or registered trademarks of their respective owners and manufacturers.